

Do you really believe it's clean?

Laboratorios
MICROKIT
★★★★★



Sight often deceives us ...




Whit KITPRO-PLUS
you can control the cleanliness of
your facilities and the health of your people.

**Colorimetric test for immediate control
of surfaces hygiene:**

- ➔ 1 year of expiry time
- ➔ Designed in easy to use and flexible sticks
- ➔ Maximum sensitivity (5 mg protein / test)
- ➔ All in a single reagent
- ➔ Immediate results
- ➔ Easy to interpret by color change
- ➔ No refrigerator need
- ➔ No apparatus are needed

COLOR CHAR

Clean Dirty Very polluted

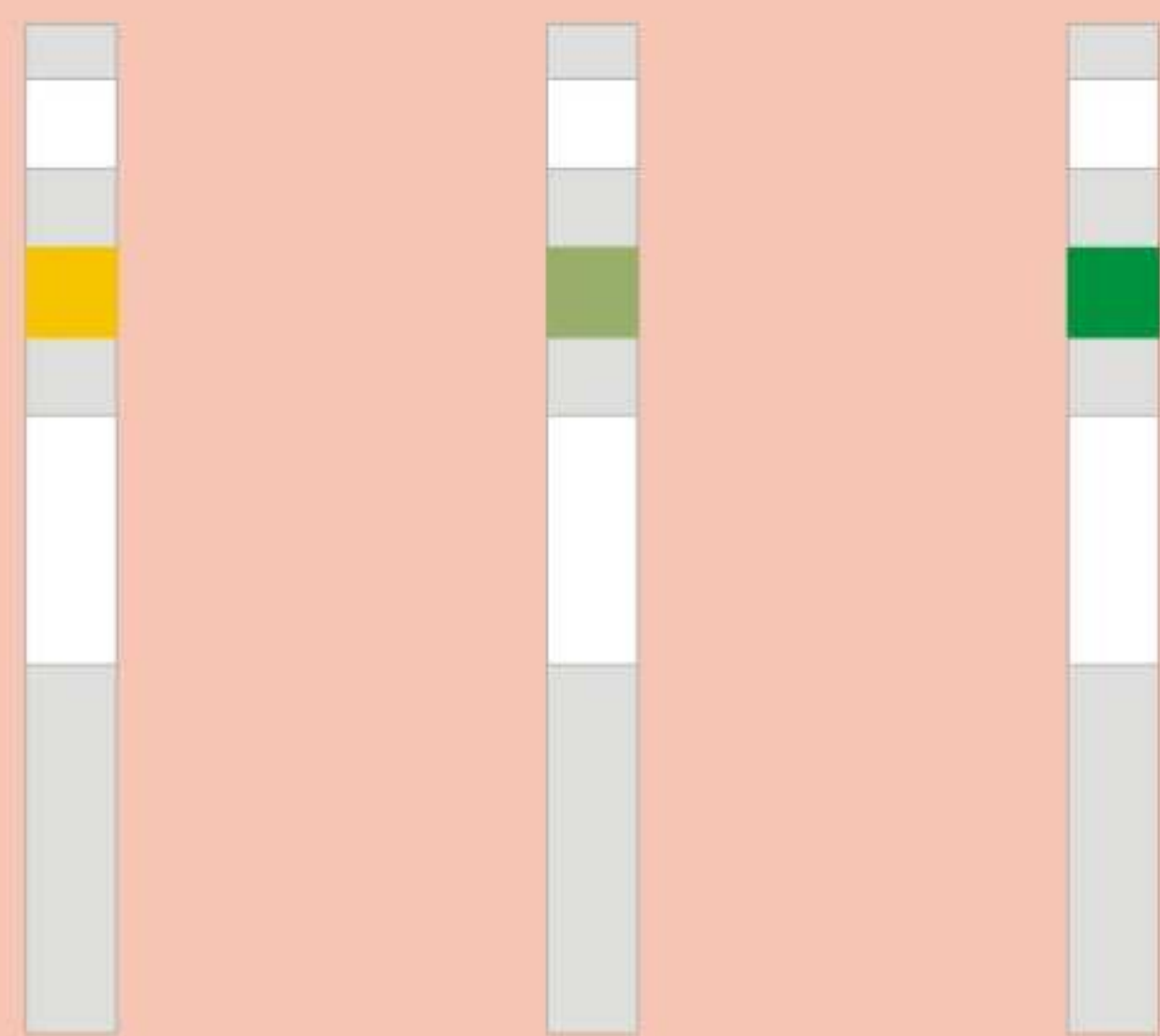

KITPRO-PLUS does not detect dust (this is already seen by your eyes). It detects the invisible dangerous dirtiness: the organic material (as proteins, such allergens). This is the more dangerous dirt, because it is the food that will multiply bacteria and fungi on your surfaces in a few hours.

KITPRO-PLUS need not qualified personnel, given the simplicity of use and interpretation.

KITPRO-PLUS instructional design allows all staff aware of the need to work with maximum hygiene.

INSTRUCTIONS FOR USE:

1. Moisten with the water spray 10 x 10 cm² of the surface. If your cleaning specifications are less stringent, do it in a 5x5, 3x3, 2x2... cm².
2. Sweep all the wetted surface with the strip, applying the finger after the yellow square, which is to be towards the surface.
3. Wait 1-60 seconds a color change from yellow to blue-green, organic and / or microbial contamination signal on the surface. The faster and stronger the color change, the greater pollution is on the surface and the faster and the greater intensity must be cleaned.

✿ Control in a simple and quick hygiene surfaces test for self-control in factories, restaurants, hotels, supermarkets, schools, hospitals, offices...

✿ Also designed to control your home hygiene, kitchens, refrigerators, tables, crockery, instrumental, showers, toilets, sinks ...